



CRYPTOCLEAR PLUS Filter Cartridges

- liquid filters
- polypropylene

CRYPTOCLEAR PLUS pleated filter cartridges have been designed specifically for the removal of *Cryptosporidium parvum* and *Giardia intestinalis* from water in the food, beverage and healthcare industries.

Extensive research, including live oocyst challenge has resulted in a graded density filtration medium that maximises loading capacity of the filters whilst accurately defining particle and oocyst retention under a variety of operating conditions.


CRYPTOCLEAR PLUS cartridges can be repeatedly sanitised using hot water, steam and a wide range of chemicals.

Features and Benefits

- Specifically designed for the reduction of *Cryptosporidium parvum* oocysts
- 0.6 and 1.0 micron retention ratings
- All polypropylene construction
- Graded density pleated media optimised dirt capacity and oocyst retention
- Independently tested with viable *Cryptosporidium parvum* oocysts



Performance Characteristics

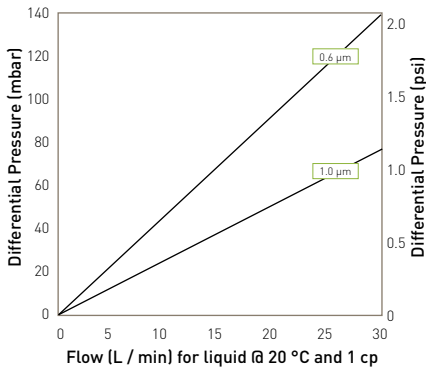

For K size for a given flow rate multiply 10" size differential pressure by 2

10" Size (250 mm) Cartridge

Specifications

Materials of Construction

- Filtration Media: Polypropylene
- Upstream Support: Polypropylene
- Downstream Support: Polypropylene
- Inner Support Core: Polypropylene
- Outer Protection Cage: Polypropylene
- End Caps: Polypropylene
- End Cap Insert: 316L Stainless Steel
- Standard o-rings/gaskets: Silicone / EPDM
- Capsule Body: Polypropylene
- Capsule Vent Seals: Silicone

Food and Biological Safety

Materials conform to the relevant requirements of 21CFR Part 177, EC1935 / 2004 and current USP Plastics Class VI - 121 °C and ISO10993 equivalents. CRYPTOCLEAR PLUS is listed in the Water Fittings and Materials Directive Part II as a WRAS Approved Product.

WRAS - Water Regulations Advisory Scheme BS6920 Test of Effect on Water Quality

Recommended Operating Conditions

Up to 70 °C (158 °F) continuous operating temperature and higher short-term temperatures during CIP to the following limits:

Temperature °C	Temperature °F	Max. Forward dP (bar)	Max. Forward dP (psi)
20	68	5.0	72.5
40	104	4.0	58.0
60	140	3.0	43.5
80	176	2.0	29.0
90	194	1.0	14.5
>100 (steam)	>212 (steam)	0.3	4.0

Capsules may be operated up to a temperature of 40 °C (104 °F) at line pressures up to 5.0 barg (72.51 psig) for liquids and 4.0 barg (58.01 psig) in air / gas.

Effective Filtration Area (EFA)

10" (250 mm) 0.57 m² (6.13 ft²)

Cleaning and Sterilisation

CRYPTOCLEAR PLUS cartridges can be repeatedly steam sterilised in situ or autoclaved at up to 142 °C (287.6 °F). They can be sanitised with hot water at up to 90 °C (194 °F) and are compatible with a wide range of chemicals. Capsules can be repeatedly autoclaved up to 135 °C (275 °F).

For detailed operational procedures and advice on cleaning and sterilisation, please contact the Technical Support Group through your usual Parker domnick hunter contact.

Retention Characteristics

The removal efficiencies of CRYPTOCLEAR PLUS cartridges have been determined from tests conducted by Thames Water Utilities Limited on live *Cryptosporidium* oocysts.

Product	Micron	Retention
CRYPTOCLEAR PLUS	0.6	>99.997%
CRYPTOCLEAR PLUS	1.0	>99.3%

Recommended Rinse Volume

Prior to use - 10 litres per 10" (250 mm) filter cartridge.

Ordering Information


Cartridges



Code	Length (Nominal)	Code	Micron	Code	Endcap (10")
B	2.5" (65 mm)	.60	0.6 µm	B	dh DOE
A	5" (125 mm)	1.0	1.0 µm	C	BF / 226 Bayonet
K	5" (125 mm)			G	Recess / 222
1	10" (250 mm)			R	BF / 222 Bayonet
2	20" (500 mm)				
3	30" (750 mm)				
4	40" (1000 mm)				

Code	Endcap (Demi)
T	TRUESEAL
Y	Demi Stub
Z	Demi A & B Std

Capsules

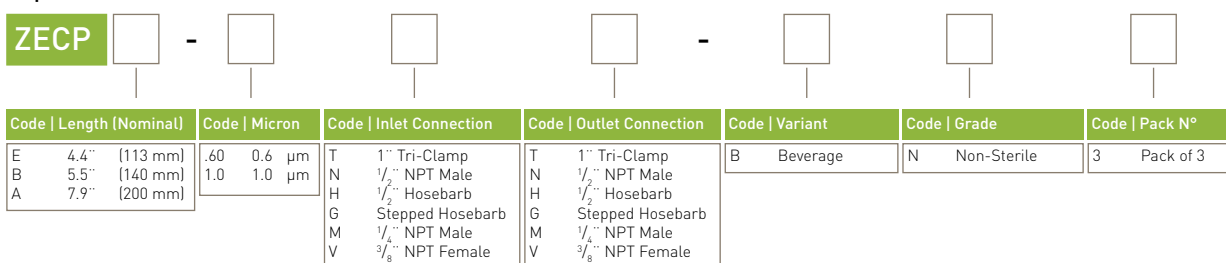

Code	Length (Nominal)	Code	Micron	Code	Inlet Connection	Code	Outlet Connection	Code	Variant	Code	Grade	Code	Pack N°
E	4.4" (113 mm)	.60	0.6 µm	T	1" Tri-Clamp	T	1" Tri-Clamp	B	Beverage	N	Non-Sterile	3	Pack of 3
B	5.5" (140 mm)	1.0	1.0 µm	N	1/2" NPT Male	N	1/2" NPT Male						
A	7.9" (200 mm)			H	1/2" Hosebarb	H	1/2" Hosebarb						
				G	Stepped Hosebarb	G	Stepped Hosebarb						
				M	1/4" NPT Male	M	1/4" NPT Male						
				V	3/8" NPT Female	V	3/8" NPT Female						

Parker domnick hunter has a continuous policy of product development and although the Company reserves the right to change specifications, it attempts to keep customers informed of any alterations. This publication is for general information only and customers are requested to contact our Process Filtration Sales Department for detailed information and advice on a products suitability for specific applications. All products are sold subject to the company's Standard conditions of sale.

12/08 Rev. 3A